


MedFrontier® FGF23 Intact


Intact Fibroblast Growth Factor 23

The better choice to measure the biologically active form of FGF23. The MedFrontier® FGF23 Intact assay kit is a reliable, reproducible, and accurate serum blood test delivering optimum results quickly.

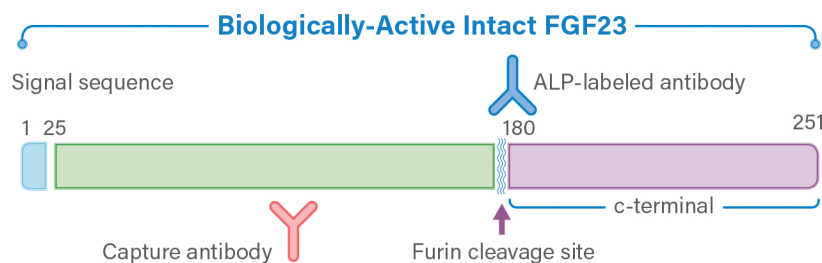


ASSAY KEY FEATURES

- ➔ Measures intact FGF23, the full-length bioactive form of FGF23
- ➔ Superior dynamic range of 10-3000 pg/mL
- ➔ Excellent reproducibility, CVs below 5%
- ➔ Microtiter plate assay that every lab can perform
- ➔ CLEIA technology
- ➔ Only 20 µL serum needed

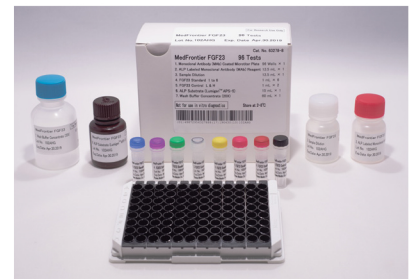

AREAS OF FGF23 RESEARCH*

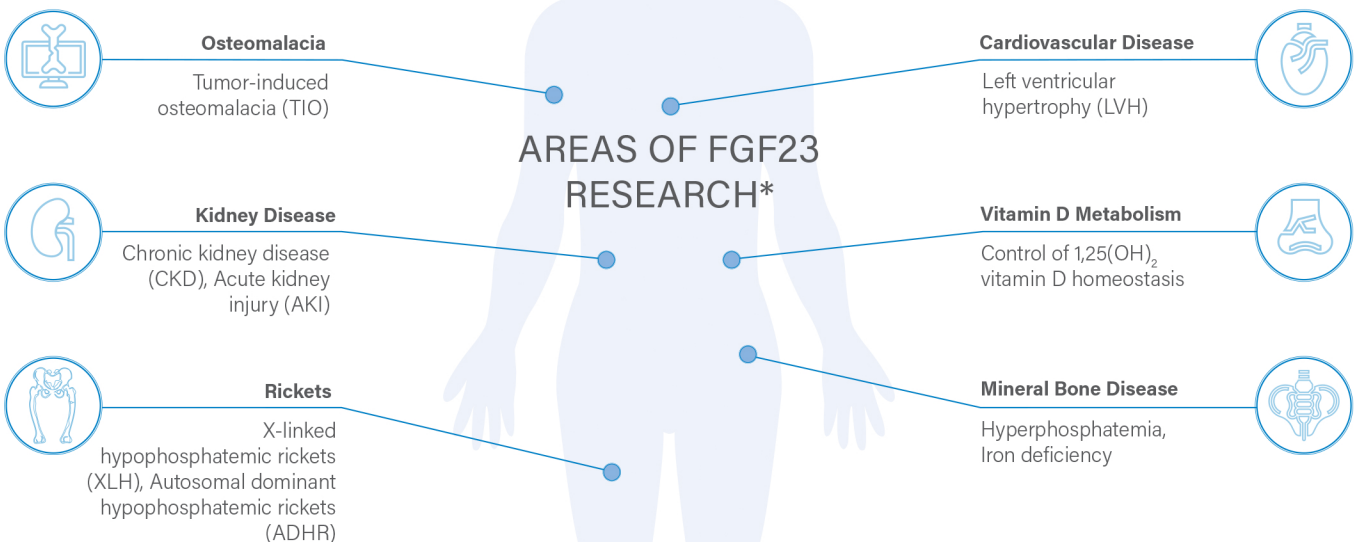

For Research Use Only. Not for use in diagnostic procedures.

*Reference: Myles Wolf et. al. Curr Opin Nephrol Hypertens. 2014 Jul; 23(4): 411-419

CHEMILUMINESCENCE ENZYME IMMUNOASSAY (CLEIA)


- ➡ CLEIA known for better sensitivity, specificity, reliability and reproducibility
- ➡ Serum intact FGF23 concentration determined using CLEIA
- ➡ Antibody-coated microtiter plate configuration for easy use
- ➡ Use of two mouse monoclonal antibodies measures only intact FGF23

ASSAY PERFORMANCE / ACCURACY & REPRODUCIBILITY

Intact FGF23 Concentration in Serum (pg/mL)	Expected Value (pg/mL)	Measured Value (pg/mL)	Accuracy (%)	Reproducibility (CV : N=10)
Low	22.8	22.7	99.6	3.1%
Medium	249.9	271.6	108.7	1.4%
High	1828.8	1946.3	106.4	2.4%

INTACT FGF23 AND VITAMIN D REGULATE PHOSPHATE HOMEOSTASIS

Intact FGF23 is the biologically active protein encoded by the FGF23 gene responsible for phosphate and Vitamin D metabolism.

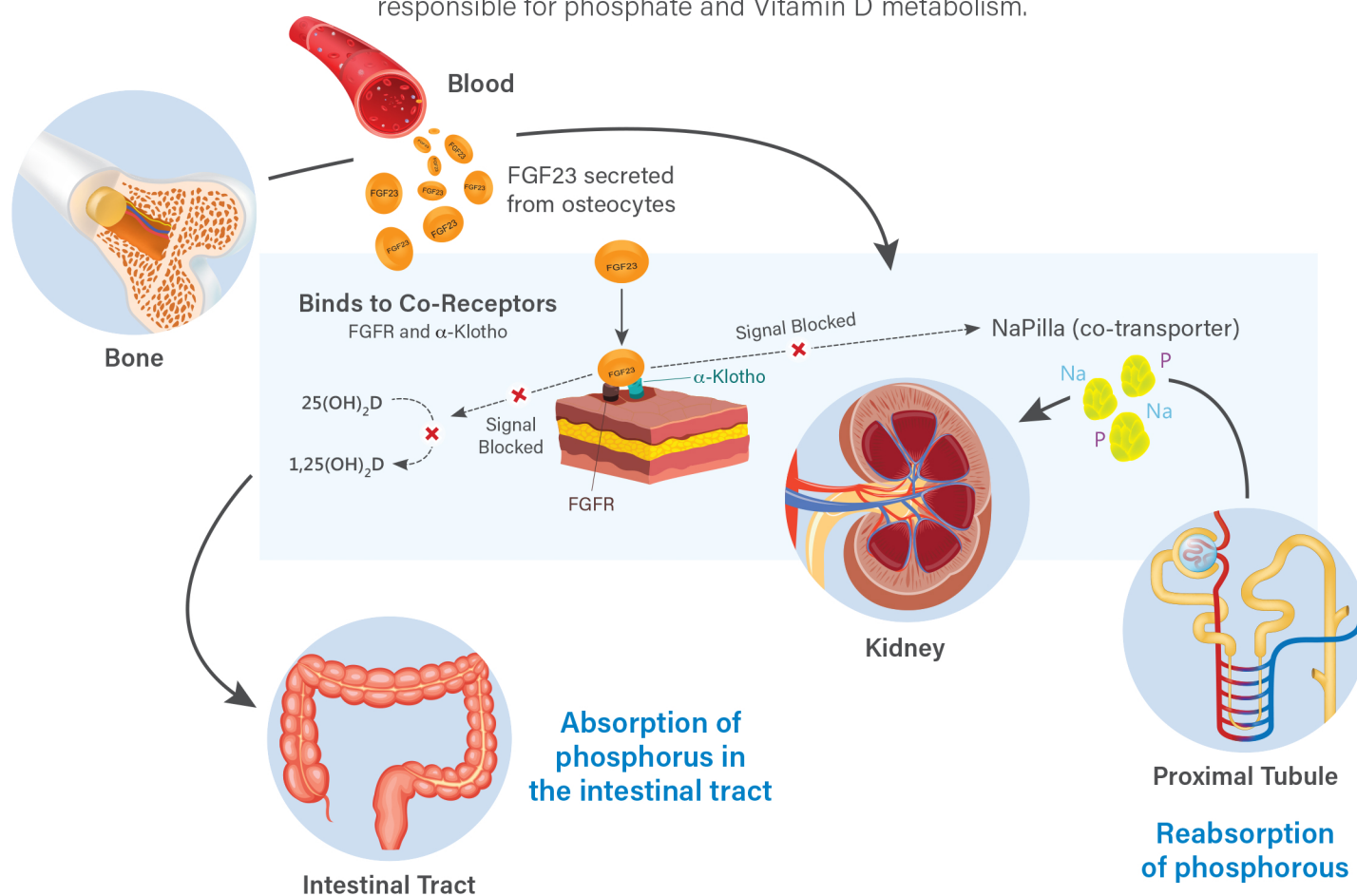

Minaris Medical Co., Ltd

1-8-10 Harumi, Chuo-ku
Tokyo 104-6004, Japan
www.minaris-medical.co.jp

For Research Use Only. Not for use in diagnostic procedures.
©2021 by Minaris Medical Co., Ltd. All rights reserved.

Eagle Biosciences

US Distributor
617-419-2019 | eaglebio.com
Kit Available Now