



EAGLE
BIOSCIENCES

High Sensitive Mouse Insulin ELISA

Catalog Number: INH11-K01

For Research Use Only. Not for use in diagnostic procedures.

v. 2 (12 OCT 24)

EAGLE BIOSCIENCES, INC.
20A Northwest Blvd., Suite 112, Nashua, NH 03063
Phone: 617-419-2019 Fax: 617-419-1110
WWW.EAGLEBIO.COM



ASSAY PRINCIPLE

This assay is a two-site enzyme-linked immunosorbent assay (ELISA). The microtiter plate is pre-coated with a monoclonal antibody against insulin. Standards and samples are added into the wells and co-incubated with a monoclonal antibody conjugated to horseradish peroxidase (HRP) enzyme. After wash step to remove any unbound substances, 3,3',5,5'-Tetramethylbenzidine (TMB) substrate is added and color develops in proportion to the amount of insulin bound initially. The assay is stopped, and the optical density of the wells is determined using a micro-plate reader. Since the increases in absorbance are directly proportional to the amount of captured insulin, the unknown sample concentration can be interpolated from a reference curve included in each assay.

MATERIALS PROVIDED

1. Microtiter Strips (96 wells), coated with a monoclonal antibody against insulin, sealed
2. 10x Wash buffer, 30 mL
3. Assay buffer, 13 mL, ready for use
4. 100x Detection antibody solution, a monoclonal antibody against insulin conjugated to horseradish peroxidase, 0.12 mL
5. Insulin standard solutions, 0 ng/mL (5 mL), 0.2 ng/mL, 0.5 ng/mL, 1.0 ng/mL, 2.0 ng/mL, 3.5 ng/mL and 7.0 ng/mL (0.15 mL each), ready for use
6. Substrate solution, 12 mL, ready for use
7. Stop solution, 12 mL, ready for use
8. Plate cover

MATERIALS REQUIRED BUT NOT PROVIDED

1. Pipettes and pipette tips
2. 96-well plate or manual strip washer
3. Buffer and reagent reservoirs
4. Paper towels or absorbent paper
5. Plate reader capable of reading absorbency at 450 nm
6. Distilled water or deionized water
7. Horizontal micro-plate shaker capable of 600 rpm

PREPARATION OF COMPONENTS

Bring all reagents and materials to room temperature before assay.

1. 1x Wash Buffer

Prepare 1x Wash buffer by mixing the 10x Wash buffer (30 mL) with 270 mL of distilled water or deionized water. If precipitates are observed in the 10x Wash buffer bottle, warm the bottle in a 37° C water bath until the precipitates disappear. The 1x Wash buffer may be stored at 2-8°C for up to one month.

2. 1x Detection Antibody

Prepare 1x Detection antibody solution by dilution of the 100x Detection antibody solution in Assay buffer, mix well. 100 µL of the 1x Detection antibody solution is required per well. Prepare only as much 1x Detection antibody solution as needed. Return the 100x Detection antibody solution to 2-8°C immediately after the necessary volume is removed.



STORAGE

The kit should be stored at 2-8°C upon receipt. Remove any unused antibody-coated strips from the micro-plate, return them to the foil pouch and re-seal. Once opened, the strips may be stored at 2-8°C for up to one month.

CALCULATION


1. Subtract the absorbance of the blank from that of standards and samples.
2. Generate a standard curve by plotting the absorbance obtained (y-axis) against insulin concentrations (x-axis). The best fit line can be generated with any curve fitting software by regression analysis. Log-log curve fitting or curve of 4- parameter can be used for calculation.
3. Determine insulin concentration of samples from standard curve.

ASSAY PROCEDURE

1. Add 10 μL of standard or sample to its respective well.
2. Add 100 μL of 1x Detection antibody solution per well.
3. Seal the plate with a plate cover. Incubate at room temperature for 90 minutes, shaking the plate at 600 rpm on a horizontal micro-plate shaker.
4. Discard the content and tap the plate on a clean paper towel to remove residual solution in each well. Add 300 μL of 1 x Wash buffer to each well. Incubate at room temperature for 20 seconds. Discard the 1x Wash buffer and tap the plate on a clean paper towel to remove residual wash buffer. Repeat the wash step for a total 4 washes.
5. Add 100 μL of Substrate solution to each well, incubate at room temperature for 15 minutes. Protect from light
6. Add 100 μL of Stop solution to each well, gently tap the plate frame for a few seconds to ensure thorough mixing.
7. Measure absorbance of each well at 450 nm immediately.

TYPICAL STANDARD CURVE

The following standard curve is provided for demonstration only. A standard curve should be generated for each assay

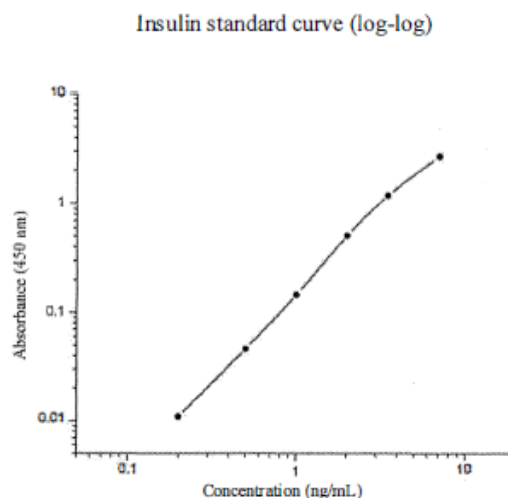



Insulin (ng/mL)	Absorbance (450 nm)	Blanked Absorbance
0	0.052	0
0.2	0.063	0.011
0.5	0.098	0.046
1.0	0.197	0.145
2.0	0.558	0.506
3.5	1.23	1.178
7	2.731	2.679

ASSAY CHARACTERISTICS

Sensitivity - The lowest insulin level that can be measured by this assay is 0.2 ng/mL.

Precision - Intra-assay Precision (Precision within an assay) C.V. <6.1 %.


Inter-assay Precision (Precision between assays) C.V. <2.9%.

Recovery - The recovery of the assay was determined by adding various amounts insulin to a sample. The measured concentration of the spiked sample in the assay was compared to the expected concentration. The average recovery was 91%.

Specificity – Percent of cross reactivity

1. Human Insulin: 100%
2. Rat Insulin: 100%

SUMMARY OF ASSAY PROCEDURE

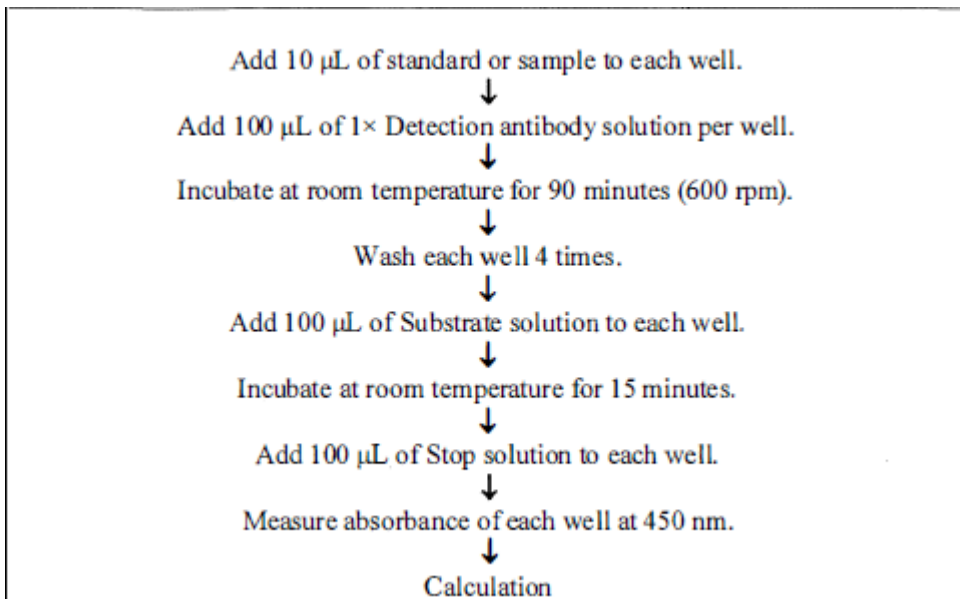



WARRANTY INFORMATION

Eagle Biosciences, Inc. warrants its Product(s) to operate or perform substantially in conformance with its specifications, as set forth in the accompanying package insert. This warranty is expressly limited to the refund of the price of any defective Product or the replacement of any defective Product with new Product. This warranty applies only when the Buyer gives written notice to the Eagle Biosciences within the expiration period of the Product(s) by the Buyer. In addition, Eagle Biosciences has no obligation to replace Product(s) as result of a) Buyer negligence, fault, or misuse, b) improper use, c) improper storage and handling, d) intentional damage, or e) event of force majeure, acts of God, or accident.

Eagle Biosciences makes no warranties, either expressed or implied, except as provided herein, including without limitation thereof, warranties as to marketability, merchantability, fitness for a particular purpose or use, or non-infringement of any intellectual property rights. In no event shall the company be liable for any indirect, incidental, or consequential damages of any nature, or losses or expenses resulting from any defective product or the use of any product. Product(s) may not be resold, modified, or altered for resale without prior written approval from Eagle Biosciences, Inc.

For further information about this kit, its application or the procedures in this kit, please contact the Technical Service Team at Eagle Biosciences, Inc. at info@eaglebio.com or at 866-411-8023.