



EAGLE
BIOSCIENCES

Mouse IFN γ Gamma ELISpot Assay

Catalog Number: RFN11-ES05

For Research Use Only. Not for use in diagnostic procedures.

v. 2 (09 SEP 24)

EAGLE BIOSCIENCES, INC.

20A Northwest Blvd., Suite 112, Nashua, NH 03063

Phone: 617-419-2019 Fax: 617-419-1110

WWW.EAGLEBIO.COM



INTENDED USE

The Eagle Biosciences ELISpot is a highly specific immunoassay for the analysis of cytokine and other soluble molecule production and secretion from T-cells at a single cell level in conditions closely comparable to the in-vivo environment with minimal cell manipulation. This technique is designed to determine the frequency of cytokine producing cells under a given stimulation and the comparison of such frequency against a specific treatment or pathological state. The ELISpot assay constitutes an ideal tool in the investigation of Th1 / Th2 responses, vaccine development, viral infection monitoring and treatment, cancerology, infectious disease, autoimmune diseases and transplantation.

Utilizing sandwich immuno-enzyme technology, Eagle Biosciences ELISpot assays can detect both secreted cytokines and single cells that simultaneously produce multiple cytokines. Cell secreted cytokines or soluble molecules are captured by coated antibodies avoiding diffusion in supernatant, protease degradation or binding on soluble membrane receptors. After cell removal, the captured cytokines are revealed by tracer antibodies and appropriate conjugates.

For further information about this kit, its application, or the procedures in this insert, please contact the Technical Service Team at Eagle Biosciences, Inc at www.EagleBio.com or at 866-411-8023.

SUMMARY

Different populations of T-cells secrete differing patterns of cytokines that ultimately lead to different immune responses. IFN γ production is a key function of Th1, CD8+ CTLs and also NK cells. IFN γ is a cytokine critical for cell mediated immunity against viral and intracellular bacterial infections and is involved in the inflammatory response following secretion via macrophage activation and stimulation of antibody secretion. IFN γ is the hallmark effector cytokine of Th1 and therefore is an excellent marker for identifying a host response to intracellular pathogens.

IFN γ is produced during infection by T cells of the cytotoxic/suppressor phenotype (CD8) and by a subtype of helper T cells, the Th1 cells. Th1 cells secrete IL-2, IL-3, TNF α and IFN γ , whereas Th2 cells mainly produce IL-3, IL-4, IL-5, and IL-10, but little or no IFN γ . IFN γ preferentially inhibits the proliferation of Th2 but not Th1 cells, indicating that the presence of IFN γ during an immune response will result in the preferential proliferation of Th1 cells.

In addition, IFN γ has several properties related to immunoregulation. IFN γ is a potent activator of mononuclear phagocytes and activates macrophages to kill tumor cells by releasing reactive oxygen intermediates and TNF α . IFN γ induces or augments the expression of MHC antigens on macrophages, T and B cells and some tumor cell lines. On T and B cells IFN γ promotes differentiation. It enhances proliferation of activated B cells and can act synergistically with IL-2 to increase immunoglobulin light-chain synthesis.

The role of IFN γ as a disease marker has been demonstrated for a number of different pathological situations including, viral infection, Autoimmune disease, transplant rejection, Diabetes and allergy.

ASSAY PRINCIPLE



A capture antibody highly specific for the analyte of interest is coated to the wells of a PVDF bottomed 96 well microtitre plate either during kit manufacture or in the laboratory. The plate is then blocked to minimize any non-antibody dependent unspecific binding and washed. Cell suspension and stimulant are added and the plate incubated allowing the specific antibodies to bind any analytes produced. Cells are then removed by washing prior to the addition of Biotinylated detection antibodies which bind to the previously captured analyte. Enzyme conjugated streptavidin is then added binding to the detection antibodies. Following incubation and washing substrate is then applied to the wells resulting in colored spots which can be quantified using appropriate analysis software or manually using a microscope.

1. 96-PVDF bottomed-well plates are first treated with 35% ethanol and then coated with capture antibody.




2. Cells are incubated in the presence of the stimulating agent. Upon stimulation they release cytokines which bind to the capture antibodies.





3. Cell removal by washing. Incubation with biotinylated detection antibody.





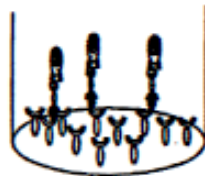

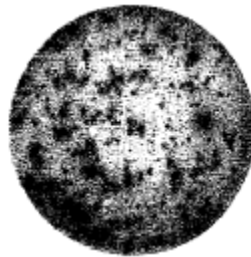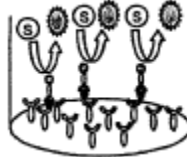

4. Any excess unbound detection antibodies is removed by washing. Incubation with streptavidin - alkaline phosphatase conjugate.



5. Any excess unbound Strep-AP is removed by washing. Incubation with BCIP/NBT. Finally BCIP/NBT reduction by alkaline phosphatase give a precipitated product which give blue/purple spots. One spot correspond to one single producing cell.



-  Capture antibody
-  Cytokines
-  Cell
-  Biotinylated detection antibody
-  Streptavidin - alkaline phosphatase conjugated
-  Substrate product



MATERIALS PROVIDED

Reagents	.001	.002	.005	Reconstitution
Pre-coated PVDF plates	1	2	5	Rehydrate with 100 μ l of PBS 1 X
Biotinylated Detection Antibody	1 (100 μ L)	1 (200 μ L)	1	Reconstitute with 0.55 ml of distilled water. Dilute prior to use
Streptavidin-Alkaline Phosphatase Conjugate	1 (10 μ L)	1 (20 μ L)	1 (50 μ L)	Dilute prior to use
Bovine Serum Albumin (BSA) – 2g	1	1	1	Dissolve to prepare dilution buffer
Ready to use BCIP/NBT – (Substrate Buffer)	1 (11ml)	1 (25ml)	2 (25ml)	Ready to use

MATERIALS PROVIDED BUT NOT PROVIDED

- Miscellaneous laboratory plastic and/or glass, if possible sterile
- Cell culture reagents (e.g. RPMI-1640, L-glutamine, FCS)
- Cell stimulation reagents (e.g. PMA, Ionomycin, Concanavalin A, mll-2)
- CO2 incubator
- Tween 20
- Phosphate Buffered Saline (PBS)

STORAGE

Store kit reagents between 2 and 8°C. Immediately after use remaining reagents should be returned to cold storage (2 to 8°C). Expiry of the kit and reagents is stated on box front labels. The expiry of the kit components can only be guaranteed if the components are stored properly, and if in the case of repeated use of one component, the reagent is not contaminated by the first handling.

SAFETY AND PRECAUTIONS FOR USE

- For research use only not to be used as a diagnostic test.
- Handling of reagents, blood specimens, PBMC, human cell lines should be in accordance with local safety procedures, e.g. CDC/NIH Health manual : "Biosafety in Microbiological and Biomedical Laboratories" 1984.
- Do not eat, drink, smoke or apply cosmetics where kit reagents are used.
- Do not pipette by mouth.
- When not in use, kit components should be stored refrigerated or frozen as indicated on vials or bottles labels.
- All reagents should be warmed to room temperature before use.
- Cover or cap all reagents when not in use.
- Do not mix or interchange reagents between different lots.
- Do not use reagents beyond the expiration date of the kit.
- Use a clean disposable plastic pipette tip for each reagent, standard, or specimen addition in order to avoid cross contamination.
- Use a clean plastic container to prepare the washing solution.
- Thoroughly mix the reagents and samples before use by agitation or swirling.



- All residual washing liquid must be drained from the wells by efficient aspiration or by decantation followed by tapping the plate forcefully on absorbent paper. Never insert absorbent paper directly into the wells.
- When pipetting reagents, maintain a consistent order of addition from well-to-well. This will ensure equal incubation times for all wells.
- BCIP/NBT substrate may cause an allergic skin reaction, caution should be taken when handling this reagent, always wear gloves.
- Follow incubation times described in the assay procedure.

REAGENT PREPARATION

1. 1X Phosphate Buffered Saline (PBS)

For 1 litre of 1 OX PBS, weigh-out:

80g NaCl,
2g KH₂PO₄,
14.4g Na₂HPO₄ ; 2H₂O.

Add distilled water to 1 litre.

***Dilute the solution to 1X before use.**

Check the pH of the 1X solution and adjust to required pH : 7.4 +/- 0.1.

2. 0.05% Tween PBS Solution (Wash Buffer)

For one plate, dilute 50 µL of Tween 20 in 100 ml of PBS 1 X.

3. 1% BSA PBS Solution (Dilution Buffer)

For one plate, dissolve 0.2 g of BSA in 20 ml of PBS 1X.

4. Detection Antibody

Reconstitute the lyophilized antibody with 0.55 ml of distilled water. Gently mix the solution and wait until all the lyophilized material is back into solution. If not used within a short period of time, reconstituted Detection Antibody should be aliquoted and stored at -20°C. In these conditions the reagent is stable for at least one year. For optimal performance prepare the reconstituted antibody dilution immediately prior to use. For one plate, dilute 100 µL of antibody into 10 ml of Dilution Buffer and mix well.

To avoid nonspecific background, it is recommended to filter the working solution using a disposable syringe and a 0.2µm filter disc.

5. Streptavidin - AP conjugate

For optimal performance, prepare the Streptavidin-AP dilution immediately prior to use. It is recommended to centrifuge the vial for a few seconds to collect all the volume at the bottom. For one plate, dilute 10 µL of Streptavidin-AP conjugate into 10 ml of Dilution Buffer and mix well. Do not keep this solution for further experiments.

To avoid nonspecific background, it is recommended to filter the working solution using a disposable syringe and a 0.2µm filter disc.

6. BCIP/NBT

The reagent is ready-to-use. It should be clear to pale yellow. If precipitates occur, filter the solution using a disposable syringe and a 0.2µm filter disc.



SAMPLE AND CONTROL PREPARATION

1. Cell Stimulation

Cells can either be stimulated directly in the antibody coated wells (Direct) or, first stimulated in 24 well plates or flask, harvested, and then plated into the coated wells (Indirect). The method used is dependent on 1) the type of cell assayed 2) the expected cell frequency. When a low number of cytokine producing cells are expected it is also advised to test them with the direct method, however, when this number is particularly high it is better to use the indirect ELISpot method. All the method steps following stimulation of the cells are the same whatever the method (direct/indirect) chosen.

2. Positive Assay Control, IFN γ production

We recommend using the following polyclonal activation as a positive control in your assay. Dilute mouse splenocytes in culture medium (e.g. RPMI 1640 supplemented with 2mM L-glutamine and 10% heat inactivated fetal calf serum) containing 0.5 μ g/ml Concanavalin A and 2ng/ml mL-2. Incubate 2 days (Indirect stimulation). Take off non adherent materials and harvest adherent cells with a cell scraper. Wash cells once. Dilute cells in culture medium supplemented with 1ng/ml PMA and 500ng/ml ionomycin (Sigma, Saint Louis, MO). Distribute 1x10⁵ to 2,5x10⁵ cells per 100 μ l in required wells of an antibody coated 96-well PVDF plates and incubate for 15-20 hours in an incubator. For other stimulators incubation times may vary, depending on the frequency of cytokine producing cells, and should be optimized in each situation.

3. Negative Assay Control

Dilute mouse splenocytes in culture medium to give an appropriate cell number (same number of unstimulated cells as stimulated sample cells) per 100 μ L with no stimulation.

4. Sample

Dilute mouse splenocytes in culture medium and stimulator of interest (i.e. Sample, Vaccine, Peptide pool or infected cells) to give an appropriate cell number per 100 μ L. Optimal assay performances are observed between 1x10⁵ and 2.5x10⁵ cells per 100 μ L. Stimulators and incubation times can be varied depending on the frequency of cytokine producing cells and therefore should be optimized by the testing laboratory.

METHOD

Assay Step		Details
1.	Addition	Add 100 μ L of PBS 1X to every well
2.	Incubation	Incubate plate at room temperature (RT) for 10 min
3.	Wash	Empty the wells by flicking the plate over a sink & gently tapping on absorbent paper
4.	Addition	Add 100 μ L of sample, positive and negative controls cell suspension to appropriate wells providing the required concentration of cells and stimulant
5.	Incubation	Cover the plate and incubate at 37 $^{\circ}$ C in a CO ₂ incubator for an appropriate length of time (15-20 hours) Note: do not agitate or move the plate during this incubation
6.	Addition	Empty the wells and remove excess solution then add 100 μ L of Wash Buffer to every well
7.	Incubation	Incubate the plate at 4 $^{\circ}$ C for 10 min



8.	Wash	Empty the wells as previous and wash the plate 3x with 100 μ L of Wash Buffer
9.	Addition	Add 100 μ L of diluted detection antibody to every well
10.	Incubation	Cover the plate and incubate at RT for 1 hour 30 min
11.	Wash	Empty the wells as previous and wash the plate 3x with 100 μ L of Wash Buffer
12.	Addition	Add 100 μ L of diluted Streptavidin-AP conjugate to every well
13.	Incubation	Cover the plate and incubate at RT for 1 hour
14.	Wash	Empty the wells and wash the plate 3x with 100 μ L of Wash Buffer
15.	Wash	Peel off the plate bottom and wash both sides of the membrane 3x under running distilled I water, once washing completely remove any excess solution by repeated tapping on absorbent paper
16.	Addition	Add 100 μ L of ready-to-use BCIP/NBT buffer to every well
17.	Development	Incubate the plate for 5-15 min monitoring spot formation visually throughout the incubation period to assess sufficient color development
18.	Wash	Empty the wells and rinse both sides of the membrane 3x under running distilled water. Completely remove any excess solution by gentle repeated tapping on absorbent paper
<p>Read Spots: allow the wells to dry and then read results. The frequency of the resulting-colored spots corresponding to the cytokine producing cells can be determined using an appropriate ELISpot reader and analysis software or manually using a microscope.</p> <p>Note: spots may become sharper after overnight incubation at 4°C in the dark Plate should be stored at RT away from direct light, but please note that color may fade over prolonged periods so read results within 24 hours.</p>		

PERFORMANCE CHARACTERISTICS

Reproducibility and Linearity

Intra-assay reproducibility and linearity were evaluated by measuring the spot development following the stimulation (PMA / Ionomycin) of 5 different mouse splenocytes concentrations, 12 repetitions. The data show the mean spot number, range and CV for the five cell concentrations.

Cells / Well	n	Mean number of spots per well	Min	Max	CV%
100000	12	TNTC	TNTC	TNTC	N/A
50000 Recommended	12	684	637	773	5.8%
25000 Recommended	12	765	739	794	2.5%
12500	12	447	408	509	7.2%
6250	12	134	113	156	10.5%



BIBLIOGRAPHY

1. Mosmann, T. R., Cherwinski, H., Bond, M. W., Giedlin, M. A., and Coffman, R. L. (1986). Two types of murine helper T cell clone. Definition according to profiles of lymphokine activities and secreted proteins. *J. Immunol.* 136, 2348-2357.
2. Gajewski, T. F., and Fitch, F. W. (1993). Anti-proliferative effect of IFN- γ in immune regulation. IFN- γ inhibits the proliferation of Th2 but not Th1 murine helper T lymphocyte clones. *J. Immunol.* 140, 4245- 4252_
3. Sastre, L., Roman, J. M., Teplow, D. 8., Dreyer, W. J., Gee, C. E., Larson, R. S., Roberts, T. M., and Springer, T. A. (1986). A partial genomic DNA clone for the alpha subunit of the mouse complement receptor type 3 and cellular adhesion molecule Mac-1. *Proc. Natl. Acad. Sci. U. S. A.* 83, 5644-5648.
4. Urban, J. L., Shepard, H. M., Rothstein, J. L., Sugarman, 8. J., and Schreiber, H. (1986). Tumor necrosis factor: a potent effector molecule for tumor cell killing by activated macrophages. *Proc. Natl. Acad. Sci. U.S. A.* 83, 5233-5237.
5. Ciampolillo, A., Guastamacchia, E., Caragiulo, L., Lollino, G., De Robertis, O., Lattanzi, V., and Giorgino, R.(1993). In vitro secretion of interleukin-1 beta and interferon-gamma by peripheral blood lymphomononuclear cells in diabetic patients. *Diabetes Res. Clin. Pract.* 21, 87-93.
6. Le thi Bich Thuy, Queen, C., and Fauci, A. S. (1986). Interferon- gamma induces light chain synthesis in interleukin 2 stimulated human B cells. *Eur. J. Immunol.* 16, 547-550.
7. Romagnani, S., Giudizi, M. G., Biagiotti, R., Almerigogna, F., Mingari, C., Maggi, E., Liang, C. M., and Moretta, L. (1986). 8 cell growth factor activity of interferon-gamma. Recombinant human interferon-gamma promotes proliferation of anti-mu-activated human B lymphocytes. *J. Immunol.* 136, 3513-3516.
8. Cunningham, A. L., Nelson, P. A, Fathman, C. G., and Merigan, T. C. (1985). Interferon gamma production by herpes simplex virus antigen-specific T cell clones from patients with recurrent herpes labialis. *J. Gen. Virol.* 66, 249-258.
9. Olsson, T. Multiple sclerosis, cerebrospinal fluid. (1994). *Ann. Neurol.* 36 Suppl, 100-102.
10. Nast, C. C., Zuo, X. J., Prehn, J., Danovitch, G. M., Wilkinson, A., and Jordan, S. C. (1994). Gamma interferon gene expression in human renal allograft fine-needle aspirates. *Transplantation* 57,498-502.
11. Suomalainen, H., Soppi, E., Laine, S., and Isolauri, E. (1993). Immunologic disturbances in cow's milk allergy, Evidence for defective interferon-gamma generation. *Pediatr. Allergy Immunol.* 4, 203-207.



Murine IFN γ ELISpot References

Huang, J. et al., Vaccine. 2006 Jun 5; 24(23): 4905-4913.
Joe, P.T. et al., J Transl Med. 2019 Jul 25;17(1):242.
Musson, J.A. et al., Infect.Immun., 2010 ;IAI.00195-10.
Patil, H.P. et al.,Vaccines (Basel). 2017 Jul 27;5(3). pii: E19.
Pollard, C. et al., Mol Ther ..,2013; 21(1): 251-9
Ramakrishna L. et al., J.Virol., 2004;78(17):9174-9189
Someya K. et al. J. Virol., 2004; 78(18):9842-9853
Tsunoda, I. et al., J Virol., 2005 ;79(23) :14640-6
Van Hoecke Let al., Nat Commun. 2018 Aug 24;9(1):3417
Van Hoecke, L. et al., Mol Ther Nucleic Acids. 2020 Dec 4; 22: 373-381. Van Hoecke, L. et al., Oncoimmunology. 2020; 9(1): 1802968.
Yan, L. et al., Infect.Immun.,2008 ;IAI.00245-08

WARRANTY INFORMATION

Eagle Biosciences, Inc. warrants its Product(s) to operate or perform substantially in conformance with its specifications, as set forth in the accompanying package insert. This warranty is expressly limited to the refund of the price of any defective Product or the replacement of any defective Product with new Product. This warranty applies only when the Buyer gives written notice to the Eagle Biosciences within the expiration period of the Product(s) by the Buyer. In addition, Eagle Biosciences has no obligation to replace Product(s) as result of a) Buyer negligence, fault, or misuse, b) improper use, c) improper storage and handling, d) intentional damage, or e) event of force majeure, acts of God, or accident.

Eagle Biosciences makes no warranties, either expressed or implied, except as provided herein, including without limitation thereof, warranties as to marketability, merchantability, fitness for a particular purpose or use, or non-infringement of any intellectual property rights. In no event shall the company be liable for any indirect, incidental, or consequential damages of any nature, or losses or expenses resulting from any defective product or the use of any product. Product(s) may not be resold, modified, or altered for resale without prior written approval from Eagle Biosciences, Inc.

For further information about this kit, its application or the procedures in this kit, please contact the Technical Service Team at Eagle Biosciences, Inc. at info@eaglebio.com or at 866-411-8023.